

METHOD 7950

ZINC (ATOMIC ABSORPTION, DIRECT ASPIRATION)

1.0 SCOPE AND APPLICATION

1.1 See Section 1.0 of Method 7000.

2.0 SUMMARY OF METHOD

2.1 See Section 2.0 of Method 7000.

3.0 INTERFERENCES

3.1 See Section 3.0 of Method 7000 if interferences are suspected.

3.2 High levels of silicon, copper, or phosphate may interfere. Addition of strontium (1,500 mg/L) removes the copper and phosphate interference.

3.3 Zinc is a universal contaminant, and great care should be taken to avoid contamination.

4.0 APPARATUS AND MATERIALS

4.1 For basic apparatus, see Section 4.0 of Method 7000.

4.2 Instrument parameters (general):

4.2.1 Zinc hollow cathode lamp.

4.2.2 Wavelength: 213.9 nm.

4.2.3 Fuel: Acetylene.

4.2.4 Oxidant: Air.

4.2.5 Type of flame: Oxidizing (fuel lean).

4.2.6 Background correction: Required.

5.0 REAGENTS

5.1 See Section 5.0 of Method 7000.

5.2 Preparation of standards:

5.2.1 **Stock solution:** Dissolve 1.000 g zinc metal (analytical reagent grade) in 10 mL of concentrated nitric acid and dilute to 1 liter with Type II water. Alternatively, procure a certified standard from a supplier and verify by comparison with a second standard.

5.2.2 Prepare dilutions of the stock solution to be used as calibration standards at the time of analysis. The calibration standards should be prepared using the same type of acid and at the same concentration as will result in the sample to be analyzed after processing.

6.0 SAMPLE COLLECTION, PRESERVATION, AND HANDLING

6.1 See Chapter Three, Section 3.1.3, Sample Handling and Preservation.

7.0 PROCEDURE

7.1 Sample preparation: The procedures for preparation of the sample are given in Chapter Three, Section 3.2.

7.2 See Method 7000, Paragraph 7.2, Direct Aspiration.

8.0 QUALITY CONTROL

8.1 See Section 8.0 of Method 7000.

9.0 METHOD PERFORMANCE

9.1 The performance characteristics for an aqueous sample free of interferences are:

Optimum concentration range: 0.05-1 mg/L with a wavelength of 213.9 nm.

Sensitivity: 0.02 mg/L.

Detection limit: 0.005 mg/L.

9.2 For concentrations of zinc below 0.01 mg/L, the furnace technique (Method 7951) is recommended.

9.3 Precision and accuracy data are available in Method 289.1 of Methods for Chemical Analysis of Water and Wastes.

10.0 REFERENCES

1. Methods for Chemical Analysis of Water and Wastes, EPA-600/4-82-055, December 1982, Method 289.1.

Method 7950
ZINC (ATOMIC ABSORPTION, DIRECT ASPIRATION)

